

HOW TO **LOAD** **LIVING IMAGE**[®] **FILES** IN *AURA* SOFTWARE:




STEP 1

Go to the "File" pull down menu.

STEP 2

Select "Load Living Image Click Folders".

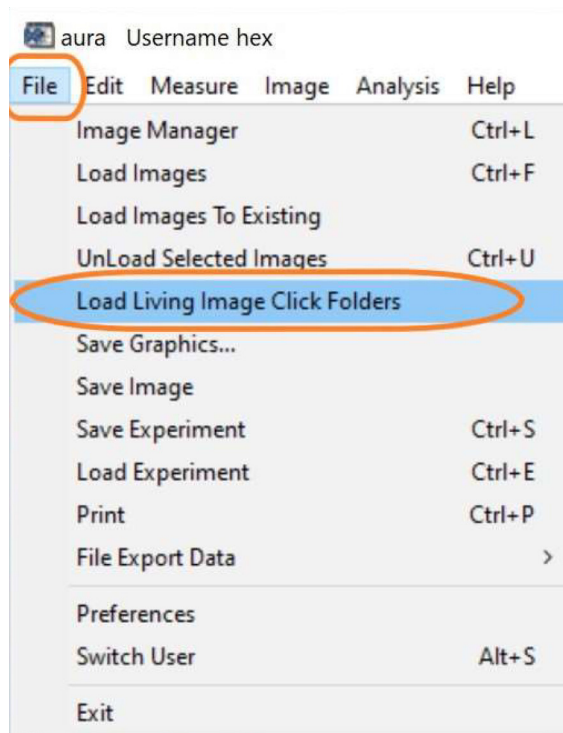


STEP 3

Locate and open the Living Image folder of interest.

Example: C:\...\Desktop\Living Image Demo Data


STEP 4A

For an individual capture, open the individual file folder. Then select the "Clickinfo.txt" file.

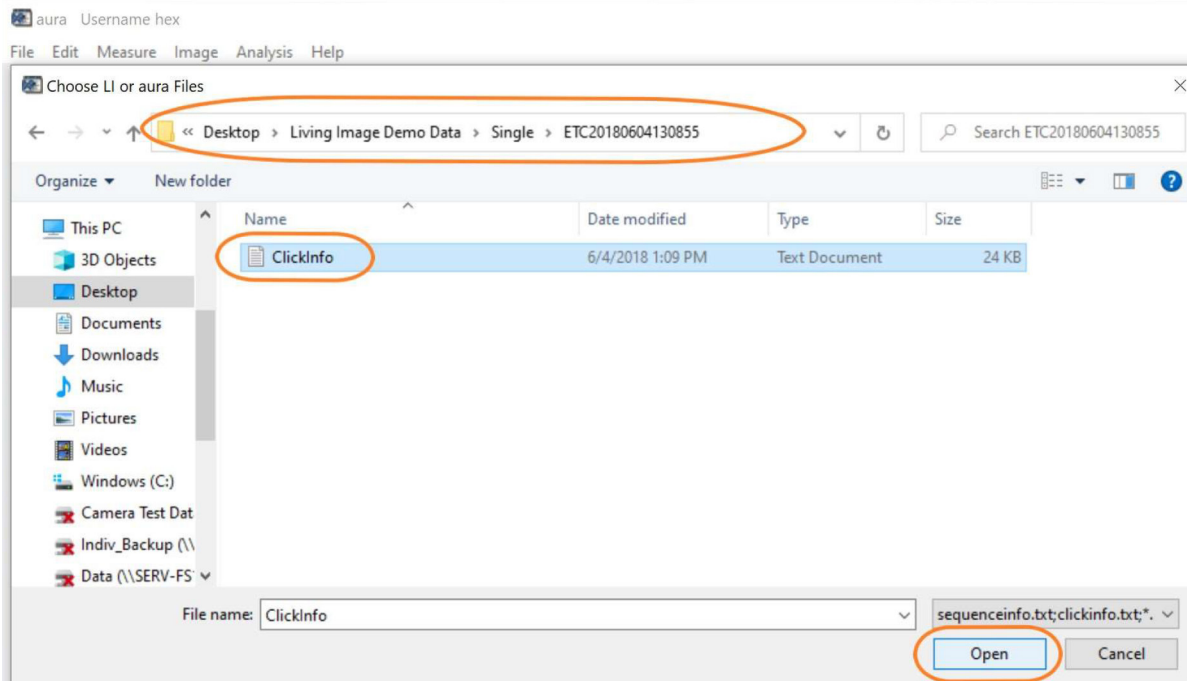

STEP 4B

For a sequence, open the sequence folder, and click on "SequenceInfo.txt".

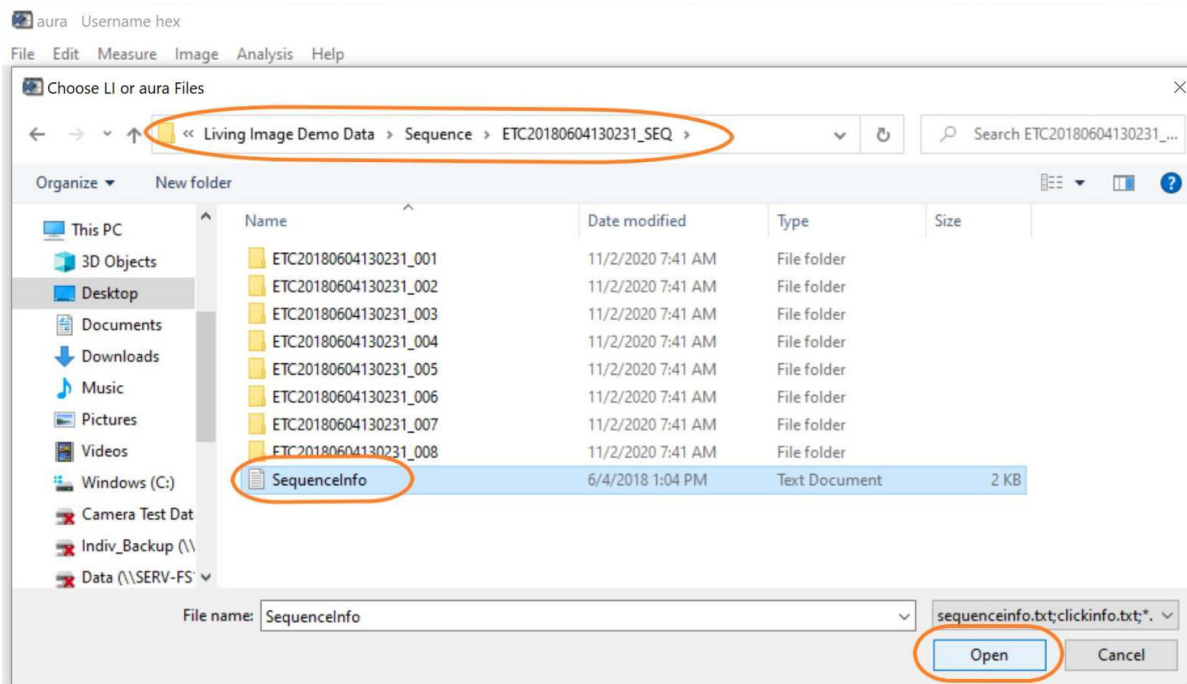

Watch Aura Analysis Tutorial Videos

